



Product Insert

Catalog # YST0019, 2X TaqMan Multiplex Real-Time qPCR Master Mix with ROX. Check our website for the various product sizes available.

TaqMan Multiplex Real-Time qPCR Master Mix is a premixed reagent solution for multiplex PCR. It has the characteristics of high amplification efficiency and high specificity. It is formulated to amplify up to 4 targets in a single reaction well and ensure excellent concordance of results in both singleplex and multiplex reactions.


Contains ROX for signal normalization.

Applications: Multiplex Real-Time qPCR.

Test Method for 20uL Reaction Volume:

Component	Volume, uL
2X TaqMan Multiplex Real-Time PCR Master Mix with ROX	10
20X TaqMan Assay (Primer/Probe Mix)	1
DNA Sample	2
Water	7
Total	20

Thermal Cycler Protocol:

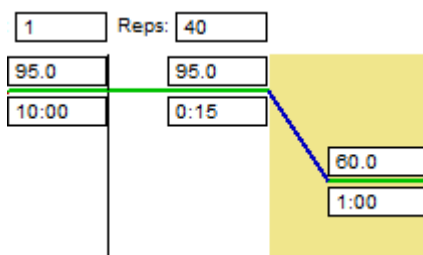

425 Boulder Court, Suite 500, Pleasanton, CA 94566, USA Phone: 1+(510) 861-2542 Email: info@sensecarebio.com

www.sensecarebio.com

FOR RESEARCH USE ONLY.