



Product Insert

YST0020, 4X TaqMan Compatible Real-Time RT-qPCR Master Mix with ROX. Check our website for the various product sizes available.

TaqMan Real-Time RT-qPCR Master Mix is a very robust 1-Step RT-qPCR master mix for RNA quantification and contains ROX for uniformity normalization.

Product Features:

- Reverse transcription and TaqMan real-time quantitative PCR are accomplished in one step for RNA quantification.
- Tolerance of inhibitors commonly found in clinical samples.

Applications: TaqMan Real-Time RT-qPCR.

Test Method for 20uL Reaction Volume:

Component	Volume, uL
2X SenseCare HotStart Taq PCR Buffer	10
20X TaqMan Assay (Primer/Probe Mix)	1
DNA Sample	2
Water	7
Total	20

Thermal Cycler Protocol:

