



Product Insert

Catalog # YST0070, 2X TaqMan Fast Versatile qPCR Master Mix with ROX. Check our website for the various product sizes available.

TaqMan Fast Versatile qPCR Master Mix with ROX has improved specificity and sensitivity for fast speed PCR instruments. It contains ROX for signal normalization. The Fast Real-Time PCR can be finished within 30-40 minutes.

Applications: Real-Time PCR for fast PCR instruments.

Test Method for 20uL Reaction Volume:

Component	Volume, uL
2X TaqMan Fast Versatile PCR Master Mix with ROX	10
20X TaqMan Assay (Primer/Probe Mix)	1
DNA Sample	2
Water	7
Total	20

Thermal Cycler Protocol:

